

Technical Release Bulletin: Morphology parameter SQA-V version 2.47.13/V-Sperm version 3.47.10

Issue date: December 30, 2007

Subject: Morphology parameters V-Sperm 3.47.10

Status/Description of failure:

- Semen Analysis Report Error: If the default setting **WHO 3rd** has been selected for morphology in the SQA-V Russian software version 2.47.13, the semen analysis report cannot be seen in the Russian V-Sperm 3.47.10. A report error will be presented.
- This error does not occur if the default setting is **WHO 4th**.

Troubleshooting to solve the problem:

Backup the data base:

1. Locate the V-Sperm folder on your computer (Normally c:\Program Files\V-Sperm).
2. Copy the entire folder to a different temporary location on your computer.

Upgrade the V-Sperm Software:


1. Close any open programs before beginning the upgrade process.
2. Insert the CD: V-Sperm Gold version 3.47.15_15 into the PC's CD-ROM. The installation process will begin automatically.
3. Follow the installation instruction (Note: The new version should be installed in the same folder it was previously installed)
4. Reboot the computer: when Installation complete: Click **OK** to restart the PC and complete the software installation

Explanation and Risks if upgrading to the new V-Sperm software version 3.47.15.15:

- The **video settings** in the new V-Sperm software version 3.47.15 are different and will need to be set-up by following the instructions below **AFTER** upgrading.

Video device that is currently installed in your system

1. After the V-sperm has been installed, turn on the V-Sperm computer and log into V-Sperm.
2. Go to Video > Video Settings and this screen will appear
3. Select these settings:
 - Video compression – DivX MPEG-4 Video Codec.
 - Video device – the video capture is currently installed.
 - Video input – Composite
 - Video size and subtype-default
 - Analog Video – pal B

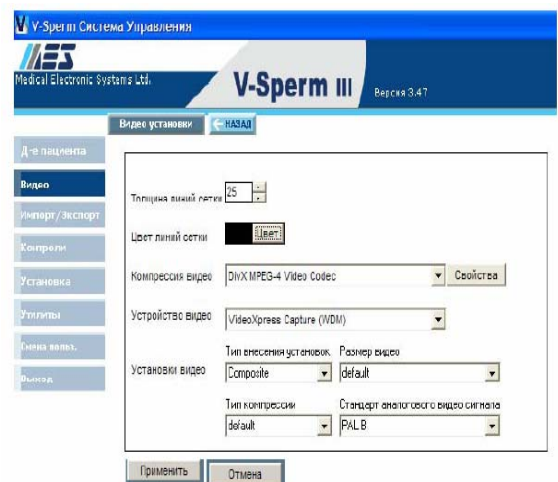

Compliance Date: Effective Immediately (30_Dec_2007)

Authority: Beni Cohen, R&D Director. beni@mes-ltd.com

Distribution: Westmedica, MES Europe



www.mes-ltd.com Tel: 972-4-637-3981 Fax: 972-4-637-3984

Technical Release Bulletin